

Moisture Analyzer

INSTRUCTION MANUAL

MS-70

MX-50

MF-50

ML-50

AND

A&D Company, Limited



Contents

1.	Safety and Compliance	3
2.	Precautions.....	6
2.1.	Installing the Analyzer.....	6
2.2.	During Use	7
2.3.	After Use and Maintaining the Analyzer	8
3.	Outline and Features.....	9
4.	Packing List and Names of each part.....	11
4.1.	Display and Keys	13
5.	Preparations	15
5.1.	Installing the Analyzer.....	15
5.2.	Setting the Clock and Calendar	16
5.2.1.	Operation	16
5.3.	Proper Operation for Precision Measurement	17
5.3.1.	Operation of the sample.....	17
5.3.2.	Operation of the analyzer	17
5.3.3.	The Glass Fiber Sheets	18
6.	Measurement Procedure	19
6.1.	Standard Mode Operation.....	19
6.1.1.	ACCURACY.....	19
6.1.2.	Operation	19
6.2.	Quick Mode Operation.....	22
6.2.1.	ACCURACY.....	22
6.2.2.	Operation	22
6.3.	Program Number	25
6.3.1.	Storing a Measurement Program to a Program Number.....	25
6.3.2.	Recalling a Measurement Program with a Program Number	25
7.	Measurement Programs	26
7.1.	List of Measurement Programs	26
7.1.1.	ACCURACY of the Standard Mode and Quick Mode	27
7.1.2.	Analyzing mode of the Automatic Mode	28
7.1.3.	Analyzing mode of the Timer Mode.....	28
7.1.4.	Drying Program (Heating Pattern and Drying Temperature).....	28
7.1.5.	Measurement Unit	29
7.2.	Procedures to Store a Measurement Program.....	30
7.2.1.	Standard Drying	30
7.2.2.	Ramp Drying	33
7.2.3.	Step Drying	37

8.	Check Function	42
8.1.	Self-Check Function	42
8.1.1.	Operation	42
8.2.	Test Sample (Sodium Tartrate Dihydrate)	43
9.	Connecting to a Printer	44
9.1.	Print samples.....	45
9.1.1.	Example To Print The Whole Data At One Time.....	45
9.1.2.	Example To Print Selected Items.....	46
9.1.3.	Explanation for Print Item	48
10.	Connecting to a Computer	50
10.1.	RS-232C Serial Interface.....	51
10.2.	Output Format	52
10.3.	Command	53
11.	Data Memory Function	54
11.1.1.	Preparation	54
11.1.2.	Output All Data at One Time	55
11.1.3.	Delete All Data at One Time	55
12.	Calibration.....	56
12.1.	Identification Number (ID No.).....	56
12.1.1.	Setting the ID Number.....	56
12.2.	Calibration of the Weighting Sensor.....	57
12.2.1.	Operation	57
12.3.	Calibration of Drying Temperature (for MS-70 and MX-50).....	59
12.3.1.	Operation	59
13.	Function Table	61
13.1.1.	Operation	62
14.	Maintenance	63
14.1.	Cleaning the Heater Unit	63
14.2.	Replacement of the Halogen Lamp	64
14.3.	Factory Settings.....	65
14.3.1.	Operation	65
14.4.	Troubleshooting.....	66
14.5.	Error Message	67
15.	Specifications.....	68
15.1.	Dimensions	69
15.2.	Accessories and Peripheral Equipment.....	70
16.	Index	71

This is a “Table of Contents preview” for quality assurance

The full manual can be purchased from our store:

<https://the-checkout-tech.com/manuals/AnD/MF-50 ML-50 MS-50 MX-70 instruction man>

And our free Online Keysheet maker:

<https://the-checkout-tech.com/Cash-Register-Keysheet-Creator/>

[HTTPS://THE-CHECKOUT-TECH.COM](https://THE-CHECKOUT-TECH.COM)