

F Series

Top Pan Balances

Operator Instructions

78102-360 Issue 2

01.04.98



Contents

- 1 Before You Begin**
 - Certificates of Conformity
- 2 The F Series**
- 3 Operating the Balance**
- 4 Balance Operating Menu**
- 5 Setting Operating Parameters**
- 6 Weighing Using Two Weighing Levels**
- 7 Interface Parameter Settings**
- 8 Utilities For Data Transfer**
- 9 Additional Functions**
- 10 Application Programs**
- 11 ISO/GLP Compliance**
- 12 Looking After Your Machine**
- 13 Appendix**

A

- Accessories, 2–4
- Additional functions, 9–1
 - blocking the keys, 9–2
 - menu access switch, 9–1
 - menu access switch on verified balances, 9–1
 - power on mode, 9–4
 - remote control, 9–3
- Adjustment, 3–3
- Animal weighing, 10–17
 - automatic start, 10–17
 - autostart storage threshold, 10–20
 - changing the number of the individual operations, 10–19
 - delayed start, 10–18
 - manual start, 10–17
- Anti-theft device, 2–7
- Application programs, 10–1
 - examples, 10–1, 10–3, 10–9, 10–12, 10–14, 10–19, 10–20
 - animal weighing/averaging, 10–17
 - counting, 10–12
 - tare memory, 10–1
 - weighing in percent, 10–6
- Averaging, 10–17

B

- Balance operating menu, 4–1
 - calling up the menu, 4–3
 - changing menu settings, 4–2
 - exiting the menu, 4–6
 - menu access switch, 4–1
 - resetting the menu, 4–7

C

- Calibration, 3–3
 - blocking calibration functions, 3–6
 - external, 3–5
 - internal, 3–4
 - sensitivity test, 3–6
- CAUTION
 - damage to your balance, 1–5
 - RFI, 1–4
- Cleaning, 12–2
- Counting, 10–12
 - changing the reference sample quantity, 10–12
 - displaying the reference sample quantity, 10–12
 - storage parameter, 10–13

D

- Data transfer utilities, 8–1
 - automatic output of application parameters, 8–3
 - automatic output of tare memory data, 8–4
 - automatic printing, 8–2
 - automatic taring (after data output), 8–3
 - data output at defined intervals, 8–3
 - data output parameter, 8–1
 - ID codes, 8–4
 - print on request, 8–3
- Description
 - connections, 2–2, 2–10
 - display, 2–14
 - general, 2–1
 - keypad, 2–16
 - options, 2–4
 - switches, 2–2, 2–10
- Display, symbols, 2–14

E

- EU verified (M) balances, 1–7
 - menu access switch, 9–1
 - specifications, 13–8
- Examples, application programs
 - animal weighing (auto start mode), 10–20
 - calculation of residual percentage weight, 10–9
 - changing the number of individual operations, 10–19
 - counting small parts, 10–14
 - net total, 10–3
 - reference sample size 20 pieces, 10–12
 - tare–net–gross weighing, 10–1
- External connections, 2–10

F

- F series, models in range, 2–2
- Function keys, 2–16

I

- Installation, 1–3
 - handling, 1–3
 - connecting peripheral devices, 1–5
 - display unit adjustment (FC161), 1–6
 - electrostatic samples, 1–5
 - legal requirements, 1–4
 - positioning, 1–4
 - radio frequency interface (RFI), 1–4
 - preparation, 2–8
- Interface parameter setting, 7–1

- ISO/GLP compliance, 11–1
 - application, 11–1
 - operating the balance, 11–2
 - printout/record (how to), 11–4
 - printout/record for adjustment procedures, 11–3
 - printout/record for application programs, 11–5

M

- Maintenance, 12–1
 - cleaning, 12–2
 - servicing, 12–1

O

- Operating requirements, 1–7
 - EU verified (M) balances, 1–7
- Operating the balance, 3–1
 - additional functions, 9–1
 - applications, 10–1
 - balance operating menu, 4–1
 - data transfer utilities, 8–1
 - interface parameters, 7–1
 - ISO/GLP compliant printouts and records, 11–1, 11–2
 - operating parameters, setting, 5–1
 - below balance weighing, 3–7
 - calibration (adjustment), 3–3
 - simple weighing, 3–2
 - taring, 3–1
 - weighing using two weighing levels, 6–1

P

- Preparing your balance, 2–8

S

- Safety, 1–1, 12–3
 - inspection, 12–3
- Servicing, 12–1
- Self test, 2–11
- Setting operating parameters, 5–1
 - adapting to ambient conditions, 5–1
 - adjustment, 5–6
 - auto zero, 5–5
 - calibration, 5–6
 - linearization, 5–6
 - stability range, 5–3
 - standard filling mode, 5–2
 - standard weighing mode, 5–2
 - tare parameter, 5–4
 - zero tracking, 5–5
- Solving problems, 2–13
- Specifications, 13–1
 - standard features, 13–1
 - verified balances, 13–8
- Switching on, 2–10
- Symbols, 1–2
 - display, 2–14

T

- Taring, 3–1
 - container taring, 3–1
 - zero the balance, 3–1

W

- Warm up time, 2–12
- WARNING
 - electric shock, 1–1
 - voltage rating, 2–10
- Weighing, 3–2
 - in percent, 10–6
 - below balance, 3–7
 - items in a container, 3–2
 - simple weighing, 3–2
 - using two weighing levels, 6–1

- Weighing in percent, 10–6
 - changing the reference percentage, 10–6
 - display parameter, 10–8
 - storage parameter, 10–7
- Weighing using two weighing levels, 6–1
 - display modes, 6–3
 - weight units, 6–1

This is a “Table of Contents preview” for quality assurance

The full manual can be purchased from our store:

[https://the-checkout-tech.com/manuals/avery berkel/FA-64 FA-113 FA-114 FA-214 FA-2](https://the-checkout-tech.com/manuals/avery%20berkel/FA-64%20FA-113%20FA-114%20FA-214%20FA-214)

And our free Online Keysheet maker:

<https://the-checkout-tech.com/Cash-Register-Keysheet-Creator/>

[HTTPS://THE-CHECKOUT-TECH.COM](https://THE-CHECKOUT-TECH.COM)