



# SERVICE MANUAL

## *Pioneer™ Balances*



# TABLE OF CONTENTS

	Page No.
<b>CHAPTER 1 GETTING STARTED</b>	
1.1 Introduction .....	1-1
1.2 Service Facilities .....	1-1
1.3 Tools and Test Equipment Required .....	1-2
1.3.1 Special Tools .....	1-2
1.3.2 Standard Tools and Test Equipment .....	1-2
1.4 Specifications .....	1-2
1.5 Balance Operation .....	1-5
1.5.1 Overview of the Controls .....	1-5
1.5.2 Overview of the Display Indicators .....	1-6
1.5.3 Power On .....	1-7
1.5.4 Power Off .....	1-7
1.5.5 Menu Setup .....	1-7
1.5.6 Menu Structure .....	1-7
1.5.7 The Load Cell .....	1-8
1.6 Connecting the RS232 Interface .....	1-9
1.7 RS232 Commands .....	1-10
<b>CHAPTER 2 TROUBLESHOOTING</b>	
2.1 Troubleshooting .....	2-1
2.1.1 General Procedures for Troubleshooting .....	2-1
2.2 Diagnostic Guide .....	2-1
2.2.1 Diagnosis .....	2-1
2.3 Error Codes .....	2-4
<b>CHAPTER 3 MAINTENANCE PROCEDURES</b>	
3.1 Preventive Maintenance .....	3-1
3.1.1 Preventive Maintenance Checklist .....	3-1
3.2 Opening the Balance .....	3-1
3.2.1 Variations in Opening Pioneer Models .....	3-2
3.3 Load Cell Removal and Disassembly .....	3-3
3.3.1 Variations in Pioneer Load Cells .....	3-3
3.3.2 Disassembly of Precision Balances with Internal Calibration .....	3-4
3.3.3 Removing Internal Calibration Weights .....	3-5
3.3.4 Servicing the Internal Calibration Mechanism .....	3-5
3.3.5 Reassembling the Internal Calibration Mechanism .....	3-6
3.3.6 Disassembly of Precision Balances with InCal Mechanism .....	3-6
3.3.7 Removing the Load Cell – Precision Balances .....	3-7
3.3.8 Load Cell Disassembly – Precision Balances .....	3-8
3.3.9 Position Sensor Board Removal/Replacement/Adjustment – Precision .....	3-13
3.3.10 Main Printed Circuit Board (PCB) Replacement – All Balances .....	3-14
3.3.11 Load Cell Reassembly – Precision Balances .....	3-15
3.3.12 Removing the InCal Mechanism from InCal Analytical Balances .....	3-19
3.3.13 Removing Internal Calibration Weights and Components .....	3-19
3.3.14 Re-Installing Internal Calibration Components .....	3-21
3.3.15 Load Cell Removal/Installation – Analytical Balances .....	3-21
3.3.16 Disassembling Analytical Load Cells .....	3-22
3.3.17 Position Sensor PC Board Removal/Replacement/Adjustment - Analytical .....	3-28
3.3.18 Reassembling Analytical Load Cells .....	3-29
3.3.19 Block Style Load Cell Removal & Replacement .....	3-34

## TABLE OF CONTENTS

	<b>Page No.</b>
3.3.20 Block Style Load Cell Disassembly & Replacement with Load Cell Kit.....	3-35
3.3.21 Progressive Disassembly of Block Style Load Cells .....	3-35
3.3.22 Internal Calibration Testing .....	3-36
3.3.23 InCal Motor Removal & Replacement .....	3-36
3.3.24 InCal Mechanism Removal.....	3-37
3.3.25 InCal Mechanism Installation .....	3-38
3.3.26 Position Sensor Assembly Removal/Installation on MFR Load Cell .....	3-39
3.3.27 Removing the Block Style Load Cell from the Base Plate.....	3-40
3.3.28 Installing the Block Style Load Cell on the Base Plate .....	3-41
3.3.29 Removal and Installation of the Block Style Load Cell PC Board.....	3-42
3.3.30 Removing the Ratio Beam on Block Style Load Cells.....	3-43
3.3.31 Installing the Ratio Beam on Block Style Load Cells.....	3-47
3.3.32 Installing the InCal Weight Arm .....	3-49
3.3.33 Installing the Load Cell PCB-Temperature Sensor-Position Sensor .....	3-51
3.2.34 Setting the Up/Down Stop Screw on Block Style Load Cells .....	3-53
 <b>CHAPTER 4 TESTING</b>	
4.1. Testing.....	4-1
4.1.1 Test Masses Required .....	4-2
4.2 Operational Test .....	4-1
4.3 Segment Display Test.....	4-1
4.4. Performance Tests .....	4-2
4.4.1 Precision Test.....	4-2
4.4.2 Off-Center Load Test.....	4-3
4.4.3 Off-Center Load Adjustment for Precision Models .....	4-4
4.4.4 Off-Center Load Adjustment for Analytical Models.....	4-5
4.4.5 Off-Center Load Adjustment for Block Style Models .....	4-6
4.4.6 Repeatability Test.....	4-7
4.4.7 Linearity Test.....	4-9
 <b>CHAPTER 5 DRAWINGS AND PARTS LISTS</b>	
5-1 Precision Top Load Balance Housing & Internal Parts.....	5-2
5-2 Round Pan Block Style InCal Balance: Housing & Internal Parts.....	5-4
5-3 Square Pan Block Style Square Pan Balance: Housing & Internal Parts .....	5-6
5.4 Precision Draft Shield Balance with Draft Shield: Housing & Parts .....	5-8
5.5 Analytical Balance: Housing & Internal Parts .....	5-10
5.6 Precision Balance Load Cell & Parts .....	5-12
5.7 Analytical Balance Load Cell & Parts .....	5-14
5.8 Block Style InCal Load Cell & Parts.....	5-26
 <b>Appendix A STANDARD CALIBRATION</b>	
A.1 Calibration.....	A-1
A.1.1 Calibration Masses .....	A-1
A.2 Span Calibration .....	A-1
A.3 Linearity Calibration .....	A-2
 <b>Appendix B SERVICE CALIBRATION</b>	
B.1 Entering the Service Menu .....	B-1
B.2 Ramp .....	B-1
B.3 Service Span Calibration .....	B-2
B.4 Service Linear Calibration.....	B-3

## TABLE OF CONTENTS

	<b>Page No.</b>
<b>APPENDIX C SOFTWARE SERVICE TOOL INSTRUCTIONS</b>	
C.1 Restore EEPROM Data .....	C-2
C.2 Replace Load Cell .....	C-3
C.3 Install new Main Printed Circuit Board.....	C-4
C.4 Update Software in the Balance .....	C-5
C.5 Diagnostics .....	C-6
C.5.1 Incal Weight Mechanism Testing .....	C-6

### LIST OF TABLES

<b>TABLE NO.</b>	<b>TITLE</b>	<b>Page No.</b>
1-1	Specifications: External Calibration Models (includes M Models) .....	1-2
1-2	Specifications: Internal Calibration Models (C & CM) .....	1-3
1-3	Specifications: Carat Models.....	1-3
1-4	Specifications: Gold Models (External Calibration) .....	1-4
1-5	Specifications: Gold Models (Internal Calibration) .....	1-4
1-6	Admissible Ambient Conditions.....	1-5
1-7	Functions of Display Controls.....	1-5
1-8	Display Indicators.....	1-6
1-9	COM 1 Interface Pin Connections.....	1-9
1-10	RS232 Command Table.....	1-10
2-1	Diagnostic Guide.....	2-2
2-2	Error Codes.....	2-4
4-1	Test Masses Required .....	4-1
4-2	Tolerances – Precision Models .....	4-2
4-3	Tolerances – Analytical Models.....	4-2
4-4	Repeatability Worksheet .....	4-8
5-1	Precision Top Load Balance Housing & Internal Parts .....	5-3
5-2	Round Pan Block Style InCal Balance: Housing & Internal Parts .....	5-5
5-3	Square Pan Block Style Square Pan Balance: Housing & Internal Parts.....	5-7
5.4	Precision Draft Shield Balance with Draft Shield: Housing & Parts.....	5-9
5.5	Analytical Balance: Housing & Internal Parts.....	5-11
5.6	Precision Balance Load Cell & Parts.....	5-13
5.7	Analytical Balance Load Cell & Parts.....	5-15
5.8	Block Style InCal Load Cell & Parts .....	5-17

### LIST OF ILLUSTRATIONS

<b>FIGURE NO.</b>	<b>TITLE</b>	
1-1	Pioneer, Carat and Gold Series Displays.....	1-5
1-2	Pioneer Display Indicators.....	1-6
1-3	Pioneer Menu Structure .....	1-7
1-4	Load Cell Assembly.....	1-8
1-5	COM 1 Connector .....	1-9
1-6	Soldering points J1 and J2.....	1-9
3-1	Top panel lifts up and slides out.....	3-2
3-2	Groove in side panel's base liner aligns with tab on Housing .....	3-2
3-3	Precision Load Cell with Incal .....	3-3
3-4	Analytical Load Cell with Incal.....	3-3
3-5	Precision Load Cell .....	3-3
3-6	Analytical Load Cell.....	3-3
3-7	Precision Block Style Load Cell with InCal.....	3-4

## TABLE OF CONTENTS

FIGURE NO.	TITLE	Page No.
3-8	Analytical Block Style Load Cell with InCal .....	3-4
3-9	Precision Balance with InCal Mechanism .....	3-4
3-10	InCal Weight Cover removed .....	3-5
3-11	Bottom Housing with Load Cell, Weight Arms and Weights removed.....	3-5
3-12	InCal Motor removed from its Housing.....	3-6
3-13	Screws securing Weight Arm assembly .....	3-6
3-14	Proper manner to grasp Load Cell when removing Weight Arm assembly.....	3-7
3-15	Cable connects to Position Sensor Board.....	3-7
3-16	Access the 3 screws to release Precision balance Load Cell from top center .....	3-7
3-17	Service Fixture, screws and washers for 150g to 310g Precision balances .....	3-8
3-18	Service Fixture, screws and washers for 510g to 4100g Precision balances .....	3-8
3-19	Precision Load Cell top .....	3-8
3-20	Precision Load Cell bottom .....	3-8
3-21	Precision Load Cell (150g to 310g) top, with Service Fixture attached.....	3-9
3-22	Service Fixture for 510g to 4100g Precision Load Cell, attached to Hanger.....	3-9
3-23	Bent Flexure .....	3-9
3-24	Position Sensor Board.....	3-10
3-25	Contact Board with ribbon cable .....	3-10
3-26	Fine wires connected to the Contact Board, and affixed to Ratio Beam .....	3-10
3-27	Contact Board taped to Ratio Beam after removal from frame .....	3-11
3-28	Hanger .....	3-10
3-29	Two screws secure Magnet Half .....	3-11
3-30	Sideways Motion Stop Plate.....	3-11
3-31	Ratio Beam clearing stop on Up/Down Stop Screw as it is lifted out .....	3-12
3-32	Magnets must be clean .....	3-12
3-33	Position Sensor Board.....	3-13
3-34	Adjust position of Sensor Board until <i>Err8.X</i> does not appear .....	3-13
3-35	Position Sensor Board.....	3-14
3-36	Installing Ratio Beam, clearing the Up/down Stop Screw .....	3-15
3-37	Precision Load Cell (150g to 310g) top, with Service Fixture attached.....	3-15
3-38	Service Fixture for 510g to 4100g Precision Load Cell, attached to Hanger.....	3-15
3-39	Sideways Motion Stop Plate.....	3-16
3-40	Two screws secure Magnet Half .....	3-16
3-41	Contact Board .....	3-16
3-42	Solder points on Position Sensor Board.....	3-16
3-43	Precision Load Cell top .....	3-17
3-44	Precision Load Cell bottom .....	3-17
3-45	Ratio Beam Flexures and Vertical Flexure.....	3-17
3-46	Insert the 3 screws to secure Precision Load Cell in Bottom Housing .....	3-17
3-47	When Position Sensor Board shows normal weight, tighten its screws .....	3-18
3-48	Analytical Balance with InCal Mechanism.....	3-19
3-49	InCal Weight Covers removed .....	3-19
3-50	Bottom Housing with Load Cell, Weight Arms and Weights removed.....	3-20
3-51	InCal Motor removed from its assembly.....	3-20
3-52	InCal Motor Cam in correct position for reassembly .....	3-20
3-53	Cable connects to the Position Sensor Board.....	3-21
3-54	On Analytical balances, the 3 screws securing Load Cell are on bottom.....	3-21
3-55	Service Fixture, screws and shims for Analytical balances.....	3-22
3-56	Bottom Mounting Plate and Flexure Arms, Analytical balance's Load Cell.....	3-22
3-57	Top Mounting Plate and Flexure Arms, Analytical balance's Load Cell.....	3-22

---

## TABLE OF CONTENTS

---

### LIST OF ILLUSTRATIONS

FIGURE NO.	TITLE	Page No.
3-58	Screws removed from Bottom Mounting Plate inserted into the Hanger.....	3-23
3-59	Shims are inserted between the Hanger and the Load Cell.....	3-23
3-60	Analytical balance Load Cell with Top Bracket visible .....	3-23
3-61	Vertical Flexure and screws, Lock Plate, washers, and bushings, removed.....	3-23
3-62	Top Flexure Arm (Analytical).....	3-23
3-63	Bottom Flexure Arm (Analytical).....	3-23
3-64	Hanger with two screws .....	3-25
3-65	Service Fixture attached to the front of Load Cell after removing Hanger .....	3-25
3-66	The Vane is secured by two screws near the rear of the Load Cell .....	3-25
3-67	The Vane removed.....	3-25
3-68	Two solder points on PSB, and two screw holes for removing the board .....	3-26
3-69	Contact Board with 3-wire ribbon cable soldered to it .....	3-26
3-70	Four screws secure the Magnet Half to the Ratio Beam .....	3-26
3-71	The Magnet Half, when removed, should be cleaned .....	3-25
3-72	A screw connects the Contact Board to the Load Cell .....	3-26
3-73	Use a smaller screw to temporarily fasten it to the Ratio Beam .....	3-26
3-74	The Sideways Motion Stop Plate and the black Up/Down Stop screw .....	3-27
3-75	Position Sensor Board (Analytical).....	3-28
3-76	Adjust Position Sensor Board vertically until <i>Err8.X</i> does not appear.....	3-28
3-77	The Ratio Beam needs to clear the groove in the Up/Down Stop Screw.....	3-29
3-78	Service Fixture is installed prior to installing Ratio Beam Flexures.....	3-29
3-79	Sideways Motion Stop Plate and Up/Down Stop Screw .....	3-29
3-80	Elongated holes on Ratio Beam Flexures go on the bottom .....	3-29
3-81	Installing the Magnet Half.....	3-30
3-82	Four screws on Magnet Half .....	3-30
3-83	Position Sensor Board.....	3-30
3-84	3-wire cable soldered to Contact Board .....	3-30
3-85	Installing the Vane.....	3-30
3-86	Screws removed from Bottom Mounting Plate inserted into the Hanger.....	3-31
3-87	Insert shims between the Hanger and the Load Cell on both sides .....	3-31
3-88	Bottom Flexure Arm .....	3-31
3-89	Top Flexure Arm.....	3-31
3-90	Vertical Flexure .....	3-31
3-91	Vertical Flexure and screws, Lock Plate, washers, and bushings, removed.....	3-31
3-92	Analytical InCal Load Cell with Top Bracket visible.....	3-32
3-93	Analytical balance Load Cell with Top Bracket visible .....	3-32
3-94	Bottom Mounting Plate on Analytical Load Cell .....	3-32
3-95	Ribbon cable connects PSB to Main PCB.....	3-32
3-96	When Position Sensor Board shows normal weight, tighten its screws .....	3-33
3-97	Block Style Load Cell .....	3-34
3-98	Service Toolset for Block Style Load Cells. PN 476000-030 .....	3-35
3-99	InCal Motor.....	3-36
3-100	InCal Motor Removal.....	3-36
3-101	Motor Coupler .....	3-36
3-102	Motor Shaft in counter-clockwise position .....	3-37
3-103	Mounting Screws holding InCal Mechanism .....	3-37
3-104	Removing the InCal Mechanism .....	3-37
3-105	InCal Mechanism Alignment.....	3-38
3-106	Centering InCal Weight .....	3-38

---

## TABLE OF CONTENTS

---

### LIST OF ILLUSTRATIONS

FIGURE NO.	TITLE	Page No.
3-107	Removing Position Sensor .....	3-39
3-108	Ratio Beam Pin Alignment .....	3-39
3-109	Base Plate Removal.....	3-40
3-110	Load Cell Mounting Components.....	3-41
3-111	Load Cell Alignment to Base Plate.....	3-41
3-112	Block Style Load Cell PC Board Item Locations .....	3-42
3-113	Block Style Load Cell PC Board Removed from Load Cell.....	3-42
3-114	Load Cell PCB-Temperature Sensor-Position Sensor assembly .....	3-43
3-115	Position Sensor .....	3-43
3-116	Removing Temperature Sensor .....	3-43
3-117	Load Cell PCB.....	3-44
3-118	Contact Board Cover is removed by squeezing the plastic post .....	3-44
3-119	Contact Board with cover holder off .....	3-44
3-120	Centering Pins.....	3-44
3-121	Centering Pins Installed .....	3-44
3-122	Contact Board Location.....	3-45
3-123	Contact Board with cover holder off .....	3-45
3-124	Loosening Ratio Beam Nuts.....	3-45
3-125	Ratio Beam Screws and Washers removed.....	3-45
3-126	Loosening InCal Weight Arm nuts.....	3-46
3-127	Weight Arm Screws and Washers removed.....	3-46
3-128	InCal Weight Arm removal .....	3-46
3-129	InCal Weight Arm removed .....	3-46
3-130	Aluminum Sleeves removed.....	3-47
3-131	Up/Down Stop Screw .....	3-47
3-132	Centering Pins removed.....	3-47
3-133	Removing the Ratio Beam .....	3-47
3-134	Ratio Beam Removed .....	3-47
3-135	Removing the Ratio Beam .....	3-47
3-136	Ratio Beam Removed .....	3-47
3-137	Removing the Aluminum Sleeves .....	3-47
3-138	Ratio Beam Weight Removal .....	3-48
3-139	Load Cell Critical Area.....	3-48
3-140	Magnet Area.....	3-48
3-141	Ratio Beam Vane Slotted Area .....	3-49
3-142	Four Ratio Beam sleeves are larger than the four Weight Arm sleeves .....	3-49
3-143	Replacing the Ratio Beam.....	3-49
3-144	Positioning the Ratio Beam over the Magnet.....	3-49
3-145	Centering Pin goes into first hole after upward bend in Ratio Beam.....	3-49
3-146	InCal Weight Arm before insertion .....	3-50
3-147	Inserting InCal Weight Arm .....	3-50
3-148	Weight Arm must be lifted slightly to insert Centering Pin.....	3-50
3-149	Proper method for tightening nuts while holding screw head stationary .....	3-50
3-150	Centering pins in Ratio Beam.....	3-51
3-151	Tightening Ratio Beam nuts.....	3-51
3-152	Up/Down Stop Screw .....	3-51
3-153	Repositioning the Contact Board and Screw.....	3-51
3-154	Load Cell PCB, Temperature Sensor and Position Sensor.....	3-52
3-155	Groove on left of PCB fits tab on Load Cell frame .....	3-52

---

## TABLE OF CONTENTS

---

### LIST OF ILLUSTRATIONS

<b>FIGURE NO.</b>	<b>TITLE</b>	<b>Page No.</b>
3-156	Temperature Sensor extends into a hole in the Magnet .....	3-52
3-157	Position Sensor on the Magnet .....	3-53
3-158	Pin centered in PSB .....	3-53
3-159	Load Cell PCB Test Connector .....	3-54
3-160	Test Connector Detail.....	3-54
3-161	Up/Down Stop Screw .....	3-54
4-1	Segment Display .....	4-1
4-2	Mass Placement Locations for Off-Center Load Test .....	4-3
4-3	Off-Center Load Adjustment Screws on Analytical models.....	4-4
4-4	Off-Center Load Adjustment Screws on Precision models .....	4-4
4-5	Off-Center Load Adjustments.....	4-4
5-1	Precision Top Load Balance Housing & Internal Parts.....	5-2
5-2	Round Pan Block Style InCal Balance: Housing & Internal Parts.....	5-4
5-3	Square Pan Block Style Square Pan Balance: Housing & Internal Parts .....	5-6
5.4	Precision Draft Shield Balance with Draft Shield: Housing & Parts .....	5-8
5.5	Analytical Balance: Housing & Internal Parts .....	5-10
5.6	Precision Balance Load Cell & Parts .....	5-12
5.7	Analytical Balance Load Cell & Parts .....	5-14
5.8	Block Style InCal Load Cell & Parts.....	5-26
C-1	Com Port Configuration Menu.....	C-1
C-2	The Software Service Tool's Restore EEPROM Tab .....	C-2
C-3	The Software Service Tool's Replace Load Cell Tab.....	C-3
C-4	The Software Service Tool's Replace PCB Tab.....	C-4
C-5	The Software Service Tool's Download Software Tab .....	C-5
C-6	The Software Service Tool's Diagnostics Tab.....	C-6



This is a “Table of Contents preview” for quality assurance

The full manual can be purchased from our store:

[https://the-checkout-tech.com/manuals/ohaus/Pioneer service manual.html](https://the-checkout-tech.com/manuals/ohaus/Pioneer%20service%20manual.html)

And our free Online Keysheet maker:

<https://the-checkout-tech.com/Cash-Register-Keysheet-Creator/>

[HTTPS://THE-CHECKOUT-TECH.COM](https://THE-CHECKOUT-TECH.COM)